
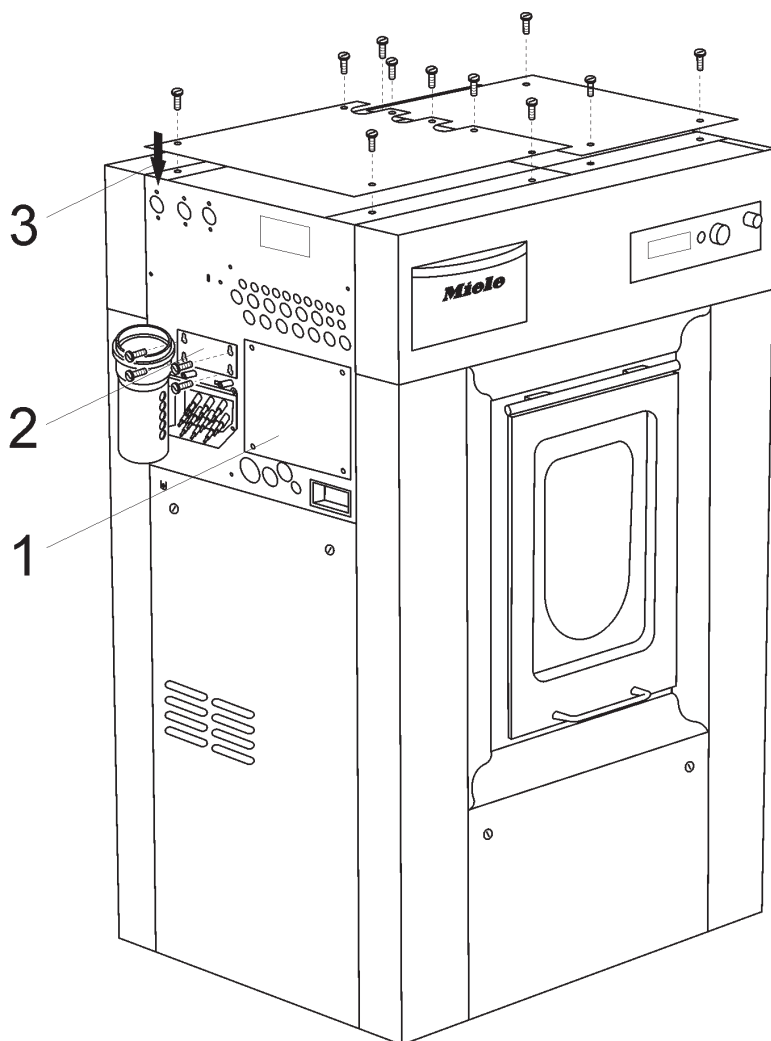
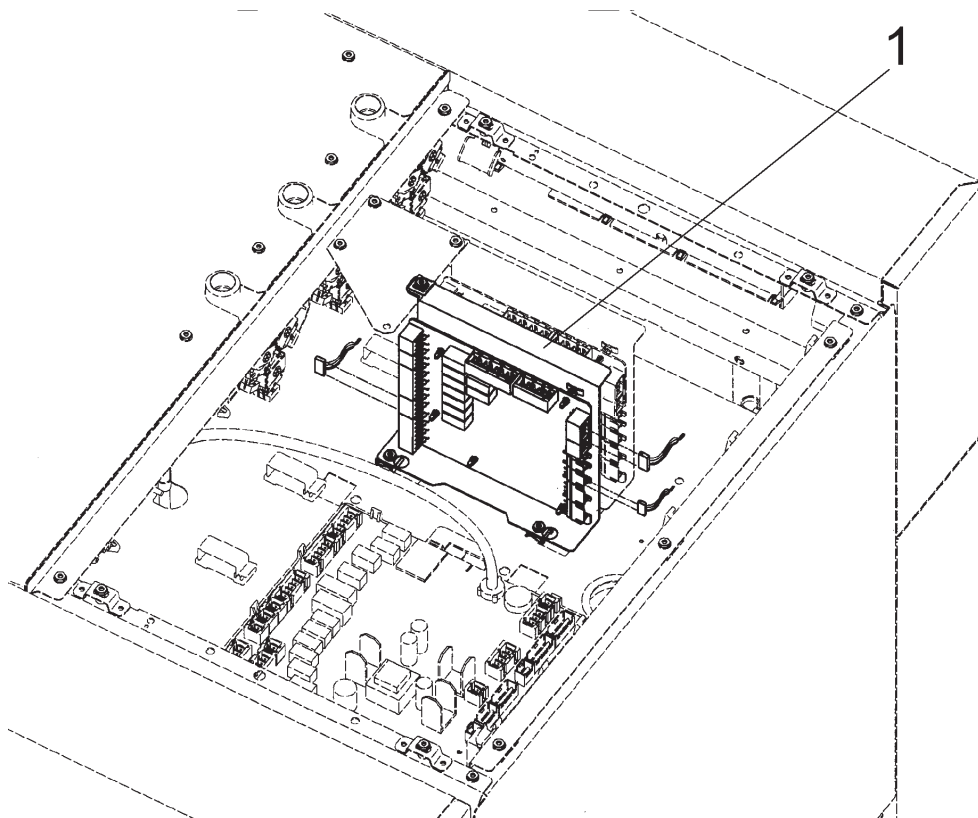
PW 6163, PW 6243, PW 6323

de	Montageanweisung Modul 1 Flüssigdosierung 1-6 Reine-Unreine-Seite	en	Fitting instructions - Module 1, liquid dispensing 1 - 6 - Clean/Soiled side	fr	Notice de montage module 1 dosage liquide 1-6 côté propre-côté contaminé
ni	Montage-instructie modul 1 vloeibarezeepdosering 1-6 schone-vuile zijde	da	Monteringsanvisning modul 1 dosering af flydende middel 1-6 ren-uren-side	sv	Monteringsanvisning: modul 1, dosering av flytande medel 1-6 ren-oren-sida
it	Istruzione di montaggio modulo 1 dosaggio liquido 1-6 lato asettico-infetto	es	Instrucciones de montaje del módulo 1 de dosificación líquida del lado descontaminado- contaminado 1-6	el	Οδηγία εγκατάστασης για μονάδα 1 δοσομέτρησης υγρού απορρυπαντικού 1-6 καθαρή-ακάθαρτη-πλευρά
am	Installation Instructions - Module 1 Liquid Dispensing 1-6 Clean-Soiled Side				


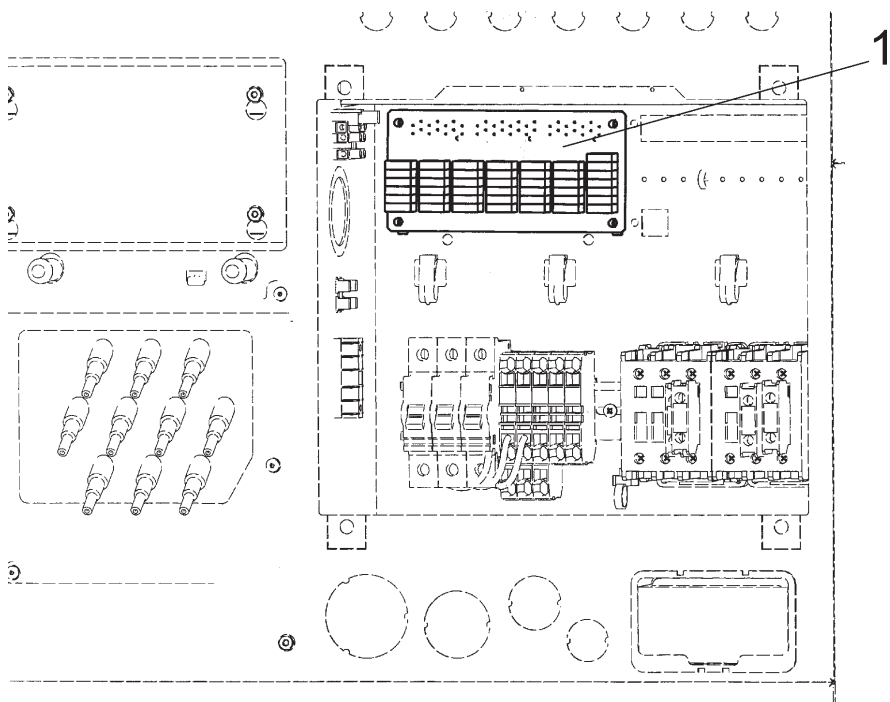
2


3

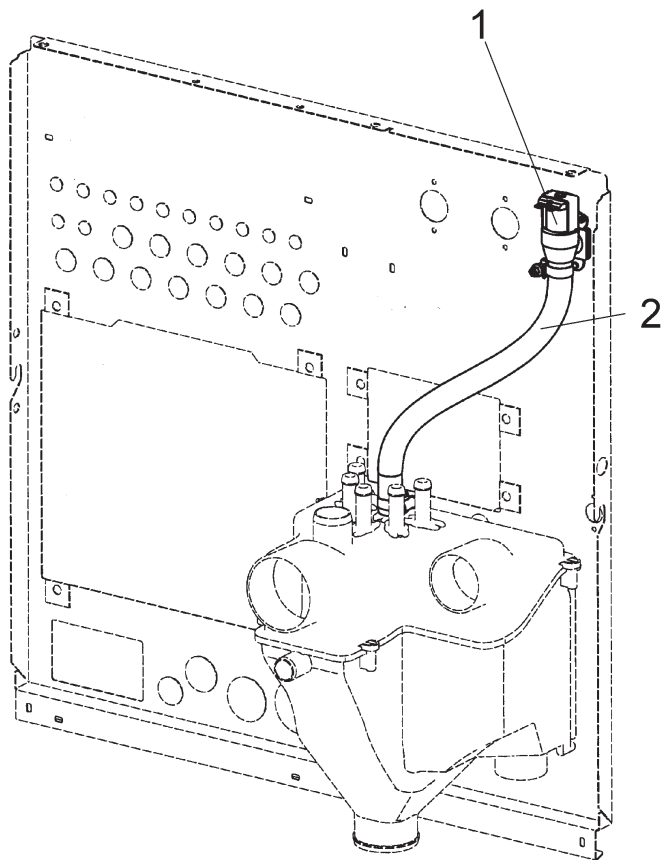4

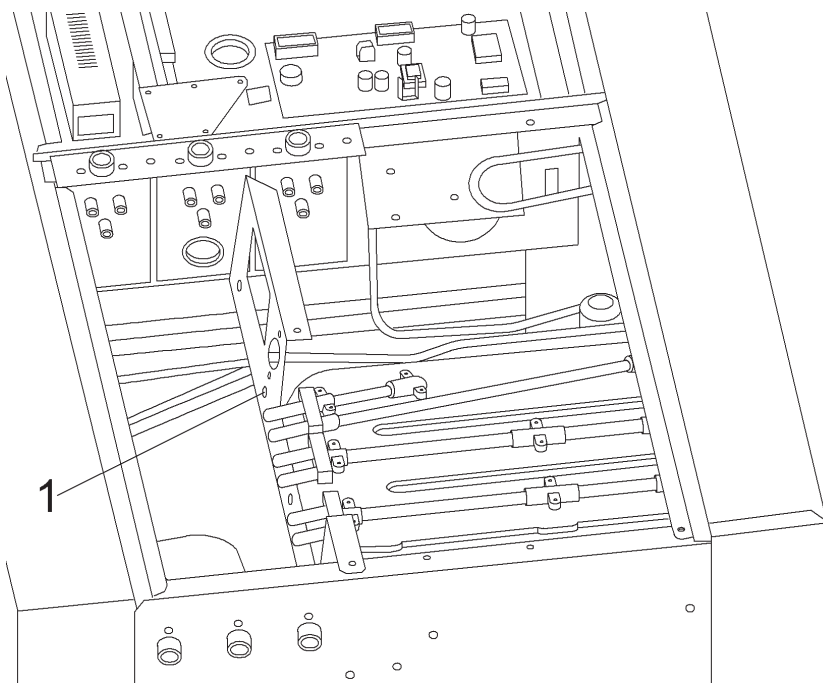5

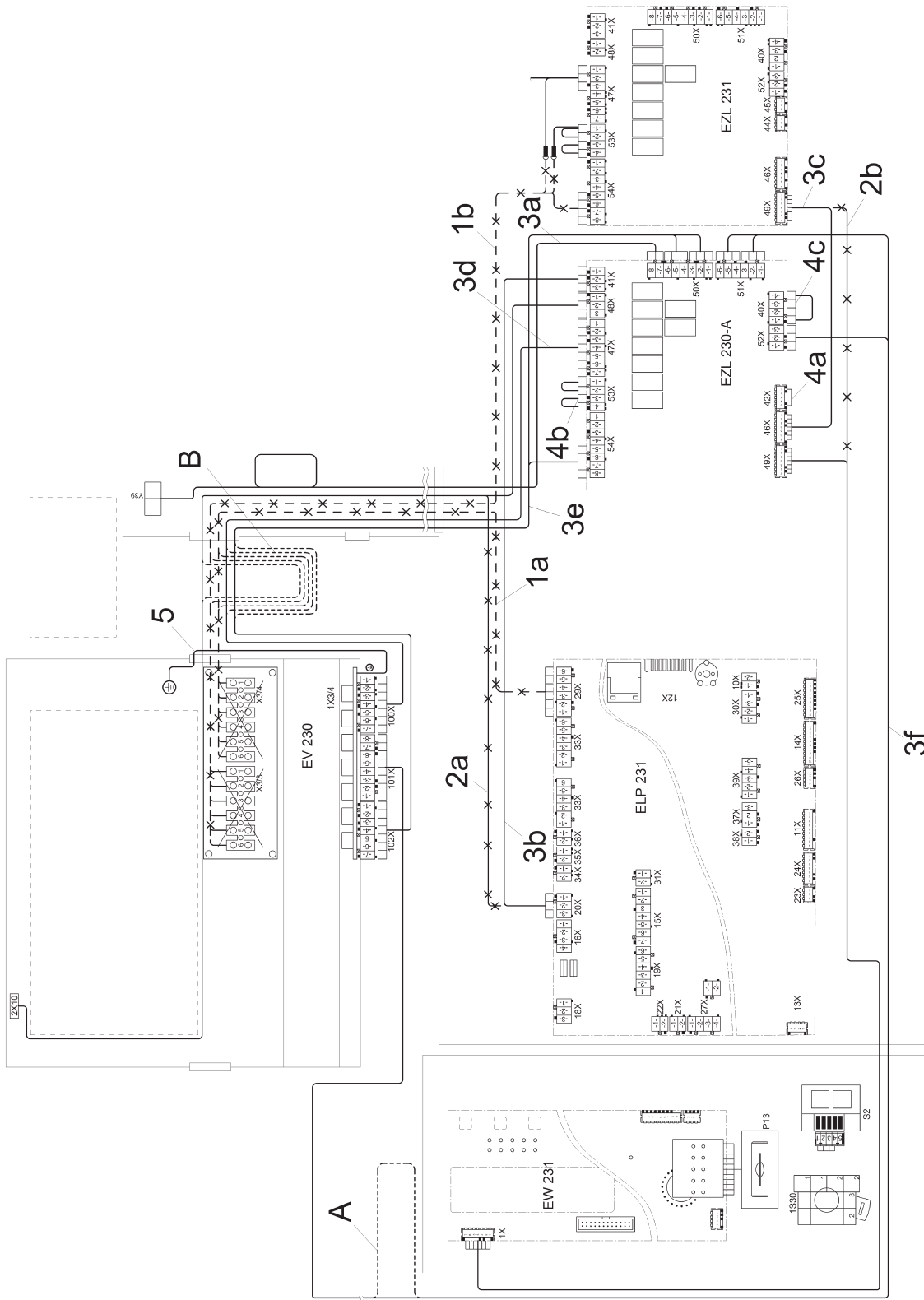de

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Note

This **service and repair work** should only be carried out by a suitably qualified electrician (with specialist training, knowledge and experience, and recent related work experience) in accordance with all appropriate local and national safety regulations.

Servicing, modification, testing and maintenance of electrical appliances should only be carried out in accordance with all appropriate legal requirements, accident prevention regulations and valid standards. All regulations of the appropriate utility supply companies and standards relating to safety (not limited to electrical safety) are to be complied with.

Danger!

Even with the machine switched off, mains voltage may be applied to some components.

Before any service work is commenced, the machine must be disconnected from the mains. Suitable measurements must be made to ensure that this is the case.

A general visual check should always be carried out.

Incorrect conversion or service work could lead to a risk of fire.

Danger!

Correct earthing function can be deactivated if casing parts are incorrectly assembled.

Correct earthing function must be ensured when refitting the casing.

Carry out appropriate electrical safety checks.

Note

The kit, Mat. no. 07285370, for PW 6163, PW 6243 and PW 6323, contains the following:

- 1 electronic module EZL 230-A, module 1 (liquid dispensing 1 - 6)
- 1 electronic module holder
- 1 coding plug MFM 1-5 (**EZL 230-A** - 42X), pin 1-5
- 1 coding plug MFM 53X (**EZL 230-A** - 53X), pin 1-2, 3-4
- 1 coding plug 40X (**EZL 230-A** - 40X), pin 1-4
- 1 adapter EV 230
- 8 spacers
- 1 wiring harness 20X-41X (**ELP 231** - 20X / **EZL 230-A** - 41X), plug housing, 3-pole, 500 mm
- 1 wiring harness 49X-46X (**EZL 231** - 49X / **EZL 230-A** - 46X), plug housing, 7-pole, 110 mm
- 1 wiring harness 47X-100X (**EZL 230-A** - 47X / **EV 230** - 100X), plug housing, 7-pole, 1820 mm
- 1 wiring harness 50X-54X-102X (**EZL 230-A** - 50X-54X / **EV 230** - 102X), plug housing, 7-4-3-pole, 1920 mm
- 1 wiring harness 51X-52X-101X (**EZL 230-A** - 51X-52X / **EV 230** - 101X), plug housing, 8-3-pole, 2600 mm
- 1 connection wire - Green-yellow 0.75 x 240

Note

The kit, Mat. no. 07285370, for PW 6163, PW 6243 and PW 6323, contains the following additional parts:

- 3 captive nuts M4
- 5 bolts M4 x 12.5
- 1 intake valve 3/4"
- 1 sticker with operating instructions information
- 1 pressure hose 330 mm
- 2 hose clips 13 - 23 mm
- 1 connection stub
- 3 cable holders 98 x 2.5
- 1 wiring harness 50X-Y39 (**EZL 230-A** - 50X / intake valve Y39), plug housing, 3-pole, 1450 mm
- 1 liquid dispensing module (1 - 6)
- 1 wiring diagram for multifunction module 1 dispensing 1 (pumps 1 - 6) and peak-load cut-out control, Mat. no. 07033893
- Fitting instructions - Module 1, liquid dispensing 1 - 6 - Clean/Soiled side, Mat. no. 07503821.

List of illustrations:

- Fig. 1: Removal of lid parts and cover plates
- Fig. 2: Fitting electronic module EZL 230-A, external view
- Fig. 3: Fitting adapter EV 230, external view
- Fig. 4: Fitting cold water connection, internal view
- Fig. 5: Cable guide connection strip
- Fig. 6: Wiring diagram EZL 230-A, module 1 (liquid dispensing 1 - 6)

- Disconnect the machine from the electrical mains and ensure power cannot be switched on again in error.
- Remove the lid parts.
- Remove the connection box cover plate (Pos. 1), liquid dispenser cover plate (Pos. 2) and intake valves cover plate (Pos. 3), see Fig. 1.

Note

If the wiring harness **ELP 231** 29X-X3/4 (external dispenser connection) is fitted, it must be removed.

If the wiring harness **EZL 231** 54X-X3/3 (peak-load cut-out connection) is fitted, it must be removed.

See Fig. 6, Pos. 1a, Pos. 1b.

- Loosen the holding plate holder.
- Slide the captive nuts M4 over the holes in the component bracket.
- Fasten the electronic module holder including the fitted additional module EZL 230-A to the component bracket with the bolts M4 x 12.5, see Fig. 2, Pos. 1.
- Tighten the holding plate holder.
- Fit the adapter EV 230 with 4 spacers on the connection box component bracket, see Fig. 3, Pos. 1.
- Connect the adapter EV 230 earth, see Fig. 6, Pos. 5.
- Fit the cold water intake valve with 2 bolts M4 x 12.5, see Fig. 4, Pos. 1.
- Fit the pressure hose 330 mm and secure it with the hose clips. Take care with the hose position on the mixer box and the position of the hose clip screws, see Fig. 4, Pos. 2.
- If there is no embossed operating instruction information under the cold water intake valve, then attach the operating instruction information sticker from the outside under the cold water intake valve.

Modify the existing plug connections as follows:

- Disconnect plug **ELP 231** - 20X and fit it in socket (module 1, liquid dispensing 1 - 6) **EZL 230-A** - 48X, see Fig. 6, Pos. 2a.
- Disconnect plug **EZL 231** - 49X and fit it in socket (module 1, liquid dispensing 1 - 6) **EZL 230-A** - 49X, see Fig. 6, Pos. 2b.

Lay and connect the wiring harnesses as follows:**Warning!**

The power and control wiring harnesses must be laid separately in order to avoid interference in the control circuit (electromagnetic compatibility).

Excess cable should be gathered such that current flows in opposite directions and tied to the cross strut in order to avoid any antenna effect, see Fig. 6, Pos. B.

- Lay the wiring harness (power) 50X-Y39 from module 1 (liquid dispensing 1 - 6) **EZL 230-A** - 50X, parallel to the existing wiring harnesses on the connection strip along the component bracket, Fig. 5, Pos. 1, to the intake valve Y39, see Fig. 6, Pos. 3a.
- Connect the cold water intake valve (Y39) in accordance with the wiring diagram.
- Lay the wiring harness (power) 20X-41X from the **ELP 231** - 20X to module 1 (liquid dispensing 1 - 6) **EZL 230-A** - 41X, see Fig. 6, Pos. 3b.
- Lay the wiring harness (control) 49X-46X from the **EZL 231** - 49X to module 1 (liquid dispensing 1 - 6) **EZL 230-A** - 46X, see Fig. 6, Pos. 3c.
- Lay the wiring harness (power) 47X-100X from module 1 (liquid dispensing 1 - 6) **EZL 230-A** - 47X, parallel to the existing wiring harnesses on the connection strip along the component bracket, Fig. 5, Pos. 1, to **EV 230** - 100X, see Fig. 6, Pos. 3d.
- Lay the wiring harness (power) 50X-54X-102X from module 1 (liquid dispensing 1 - 6) **EZL 230-A** - 50X-54X, parallel to the existing wiring harnesses on the connection strip along the component bracket, Fig. 5, Pos. 1, to **EV 230** - 102X, see Fig. 6, Pos. 3e.

Note

Lay the wiring harness (control) 51X-52X-101X from module 1 (liquid dispensing 1 - 6) **EZL 230-A** - 51X-52X, parallel to the existing wiring harnesses along the fascia panel, to **EV 230** - 101X, see Fig. 6, Pos. 3f.

Excess cable should be gathered such that current flows in opposite directions and tied in order to avoid any antenna effect, see Fig. 6, Pos. A.

- Connect the electronic module **ELZ 230-A** and adapter **EV 230** - 1X3/4 in accordance with the wiring diagram.

Check the coding plugs and modify them if necessary as follows:

- Check coding plug MFM 1-5, module 1 (liquid dispensing 1 - 6) **EZL 230-A** - 42X for “Coding 1-5”, see wiring diagram.
See Fig. 6, Pos. 4a.
- Check coding plug MFM 53X, module 1 (liquid dispensing 1 - 6) **EZL 230-A** - 53X for “Pin 1-2 and 3-4”, see wiring diagram.
See Fig. 6, Pos. 4b.
- Check coding plug 40X, module 1 (liquid dispensing 1 - 6) **EZL 230-A** - 40X for “Pin 1-4”, see wiring diagram.
See Fig. 6, Pos. 4c.

Register module 1, liquid dispensing 1 - 6 - Clean/Soiled side via the service mode.

- Service level ⇒ Main menu ⇒ Service settings ⇒ Module registration ⇒ EZL with coding 1 ⇒ Register module? ⇒ OK button ⇒ Activate disk symbol to save setting.
- Dispenser display activation.
Service level ⇒ Main menu ⇒ Service settings ⇒ Basic settings ⇒ Dispensing ⇒ Dispensing: No ⇒ Yes ⇒ OK button ⇒ Back button (to go back one level) ⇒ Activate disk symbol to save setting.

Note

If the dispensing display is activated, the operator can activate and deactivate dispensing. If the dispensing display is not activated, dispensing activation and deactivation is controlled by the programme in operation only.

- To quit the service mode, switch off the machine.

Connect the dispenser pumps in accordance with the wiring diagram and calibrate them as follows:

- If the pump capacity (flow rate) is known, enter the figure directly in the supervisor level as follow (e.g. for dispenser pump 1).
Supervisor level ⇒ Main menu ⇒ Settings ⇒ Dispensing capacity ⇒ Dispenser pump 1: No ⇒ml/min
⇒ OK button ⇒ Back button (to go back one level) ⇒ Activate disk symbol to save setting.
- If the pump capacity (flow rate) is unknown, calibrate the dispenser pump in the supervisor level as follows (e.g. for dispenser pump 1).
Supervisor level ⇒ Main menu ⇒ Settings ⇒ Dispenser calibration ⇒ Dispenser pump 1:
Place the dispenser pump suction tube in a container with a level mark.
Fill the container to the level mark.
Press button 0/1 (dispenser pump operates for 60 s).
Use a measuring vessel to refill the container to the level mark.
Note the dispensed quantity via the measuring vessel and enter this figure in the display (V:.....ml).
Activate the disk symbol to save the setting.
- Register the empty signal as follows (e.g. for dispenser pump 1).
Supervisor level ⇒ Main menu ⇒ Settings ⇒ Dispenser empty signal ⇒ Dispenser pump 1: No ⇒ Yes ⇒ OK
button ⇒ Back button (to go back one level) ⇒ Activate disk symbol to save setting.
- Check dispensing in a programme and carry out adjustments if necessary.
- Carry out a functional check.
- Carry out a data backup via the service mode.
Service level ⇒ Backup ⇒ Backup data to EW (selection module) ⇒ Overwrite existing data? ⇒ Activate
disk symbol to overwrite EW data.
- To quit the service mode, switch off the machine.

Carry out final work:

- Check the machine for leaks.
- Refit the lid parts and cover plates.
- Carry out appropriate electrical safety checks.
- Reconnect the machine to the electric mains supply.

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Note

Service and repair work should only be carried out by a suitably qualified electrician (with specialist training, knowledge and experience, and recent related work experience) in accordance with all appropriate local and national safety regulations.

Servicing, modification, testing and maintenance of electrical appliances should only be carried out in accordance with all appropriate legal requirements, accident prevention regulations and valid standards.

All regulations of the appropriate utility supply companies and standards relating to safety (not limited to electrical safety) are to be complied with.

Danger!

Even with the machine switched off, supply voltage may be applied to some components.

Before starting any service work, the machine must be disconnected from the power supply. Suitable measurements must be made to ensure that this is the case.

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Danger!

Grounding can be made ineffective if parts are reassembled incorrectly.

When reassembling the casing parts, it must be ensured that the grounding function operates correctly.

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Note

This kit, mat. no. 07285370 for PW 6163, PW 6243 and PW 6323, contains:

- 1 electronic EZL 230-A, module 1 (liquid dispensing 1-6)
- 1 electronics mounting bracket
- 1 keyed plug MFM 1-5 (**EZL 230-A** – 42X), pin 1-5
- 1 keyed plug MFM 53X (**EZL 230-A** – 53X), pin 1-2, 3-4
- 1 keyed plug 40X (**EZL 230-A** – 40X), pin 1-4
- 1 adapter EV 230
- 8 spacers
- 1 wiring harness 20X-41X (**ELP 231** – 20X / **EZL 230-A** – 41X), plug housing 3-pin, 500 mm
- 1 wiring harness 49X-46X (**EZL 231** – 49X / **EZL 230-A** – 46X), plug housing 7-pin, 110 mm
- 1 wiring harness 47X-100X (**EZL 230-A** – 47X / **EV 230** – 100X), plug housing 7-pin, 1820 mm
- 1 wiring harness 50X-54X-102X (**EZL 230-A** – 50X-54X / **EV 230** – 102X), plug housing 7-4-3 pin, 1920 mm
- 1 wiring harness 51X-52X-101X (**EZL 230-A** – 51X-52X / **EV 230** – 101X), plug housing 8-3 pin, 2600 mm
- 1 green/yellow ground wire 0.75x240

Note

This kit, mat. no. 07285370 for PW 6163, PW 6243 and PW 6323, also contains:

- 3 M4 nuts
- 5 M4x12.5 hex screws
- 1 3/4" intake valve
- 1 information sticker
- 1 pressure hose, 330 mm
- 2 hose clamps, 13-23 mm
- 1 adapter
- 3 cable ties, 98x2.5
- 1 wiring harness 50X-Y39 (**EZL 230-A** – 50X / intake valve Y39), plug housing 3-pin, 1450 mm
- 1 accessory module for liquid dispensing (1-6)
- 1 wiring diagram, multifunction module 1 dispensing 1 (pumps 1-6) and peak-load-control, mat. no. 07033893
- These instructions, "Module 1 liquid dispensing 1-6 clean-soiled side", mat. no. 07503821.

List of illustrations:

- Fig. 1, removing lid and casing
- Fig. 2, installing electronic EZL 230-A, external view
- Fig. 3, installing adapter EV 230, external view
- Fig. 4, installing cold-water connection, internal view
- Fig. 5, cable routing
- Fig. 6, schematic EZL 230-A, module 1, liquid dispensing 1-6

- Disconnect the machine from the electrical supply and ensure that power cannot be switched on again in error.
- Remove the lid from the appliance and the cover from the water side.
- Remove the covers from the connection box (Pos. 1), liquid dispensing (Pos. 2) and the intake valves (Pos. 3); see Fig. 1.

Note

Wiring harness **ELP 231** 29X-X3/4 (outside dispensing connection) is no longer necessary.

Wiring harness **EZL 231** 54X-X3/3 (peak-load connection) is no longer necessary.

See Fig. 6, Pos. 1a, Pos. 1b.

- Remove the electronics mounting bracket.
- Line up the provided M4 nuts over the holes on the chassis.
- Secure the electronics bracket, including the installed additional electronic EZL 230-A, with the M4x12.5 hex screws to the chassis; see Fig. 2, Pos. 1.
- Secure the bracket.
- Secure the EV 230 adapter with the 4 spacers to the connection box chassis; see Fig. 3, Pos. 1.
- Connect the EV 230 adapter ground connection; see Fig. 6, Pos. 5.
- Install the cold-water intake valve using 2 of the supplied M4x12.5 screws; see Fig. 4, Pos. 1.
- Install the 330 mm pressure hose using the provided hose clamps. Pay attention to the hose position on the mixing box, as well as the direction of the hose clamps' screws; see Fig. 4, Pos. 2.
- If there is no user information already imprinted underneath the cold-water intake valve, attach the provided user information sticker on the outside of the appliance, underneath the cold-water valve.

Modify plug connections:

- Disconnect plug **ELP 231** – 20X and connect it to **EZL 230-A** – 48X; see Fig. 6, Pos. 2a.
- Disconnect plug **EZL 231** – 49X and connect it to **EZL 230-A** – 49X; see Fig. 6, Pos. 2b.

Connect wiring harnesses:**Warning!**

The load- and control-current wiring harnesses are disconnected from one another in order to avoid interference signals in the control leads (electromagnetic compatibility).

Lay excess wiring in the opposite direction and secure, in order to avoid antenna-like behavior; see Fig. 6, Pos. B.

- Lay wiring harness 50X-Y39 (load current) from module 1 (liquid dispensing 1-6) **EZL 230-A** – 50X, parallel to the existing wiring harnesses alongside the chassis ground wire (Fig. 5, Pos. 1), to intake valve Y39; see Fig. 6, Pos. 3a.
- Connect the cold-water intake valve (Y39) according to the wiring diagram.
- Lay wiring harness (load current) 20X-41X from **ELP 231** – 20X to module 1 (liquid dispensing 1-6) **EZL 230-A** – 41X; see Fig. 6, Pos. 3b.
- Lay wiring harness (control current) 49X-46X from **EZL 231** – 49X to module 1 (liquid dispensing 1-6) **EZL 230-A** – 46X; see Fig. 6, Pos. 3c.
- Lay wiring harness (load current) 47X-100X from module 1 (liquid dispensing 1-6) **EZL 230-A** – 47X, parallel to the existing wiring harnesses alongside the chassis ground wire (Fig. 5, Pos. 1), to **EV 230** – 100X; see Fig. 6, Pos. 3d.
- Lay wiring harness (load current) 50X-54X-102X from module 1 (liquid dispensing 1-6) **EZL 230-A** – 50X-54X, parallel to the existing wiring harnesses alongside the chassis ground wire (Fig. 5, Pos. 1) to **EV 230** – 102X; see Fig. 6, Pos. 3e.

Note

Lay wiring harness (control current) 51X-52X-101X from module 1 (liquid dispensing 1-6) **EZL 230-A** – 51X-52X, parallel to the existing wiring harnesses to the fascia panel, to **EV 230** – 101X; see Fig. 6, Pos. 3f.

Lay excess wiring in the opposite direction and secure, in order to avoid antenna-like behavior; see Fig. 6, Pos. A.

- Connect electronic **ELZ 230-A** and adapter **EV 230** – 1X3/4 in accordance with the wiring diagram.

Check keyed plugs and swap if necessary:

- Check keyed plug MFM 1-5, module 1 (liquid dispensing 1-6) **EZL 230-A** – 42X for “key 1-5”; see the wiring diagram.
See Fig. 6, Pos. 4a.
- Check keyed plug MFM 53X, module 1 (liquid dispensing 1-6) **EZL 230-A** – 53X for “pins 1-2 and 3-4”; see the wiring diagram.
See Fig. 6, Pos. 4b.
- Check keyed plug 40X, module 1 (liquid dispensing 1-6) **EZL 230-A** – 40X for “pins 1-4”; see the wiring diagram.
See Fig. 6, Pos. 4c.

Register the liquid dispensing module:

- Service mode ⇒ Main menu ⇒ Service settings ⇒ Module registration ⇒ EZL with coding 1 ⇒ Register module? ⇒ OK button ⇒ Diskette symbol to save the registration.
- Switch on dispensing display.
Service mode ⇒ Main menu ⇒ Service settings ⇒ Basic settings ⇒ Dispensing ⇒ Dispensing: No ⇒ Yes ⇒ OK button ⇒ Back button (go back one level) ⇒ Diskette symbol to save.

Note

Through the switched-on dispensing display the user can select/deselect dispensing. When the display is switched off, the user can only view if dispensing is on for the current program and how much is dispensed.

- Switch the machine off to exit service mode.

Connect dispensing pumps as per wiring diagram and calibrate:

- If the flow rate of the dispensing pumps is known, enter this value directly in programming mode (e.g., dispensing pump 1).
Programming mode ⇒ Main menu ⇒ Settings ⇒ Dispensing flow rate ⇒ Dispensing pump 1:
No ⇒mL/min ⇒ OK button ⇒ Back button (go back one level) ⇒ Diskette symbol to save the flow rate.
- If the flow rate of the dispensing pumps is not known, calibrate the dispensing pumps in programming mode (e.g., dispensing pump 1).
Programming mode ⇒ Main menu ⇒ Settings ⇒ Calibrate dispensing ⇒ Dispensing pump 1
Insert the dispensing pump intake hose into a container with a “fill” mark.
Fill the container to the “fill” mark.
Press the 0/1 button (dispensing pump runs for 60 seconds).
Use a beaker to refill the container to the “fill” mark.
Check the dispensed amount in the beaker and record in the display (V:.....mL).
Diskette symbol to save the dispensed amount.
- Register an “empty” status message (e.g., dispensing pump 1).
Programming mode ⇒ Main menu ⇒ Settings ⇒ “Empty” message for dispensing ⇒ Dispensing pump 1:
No ⇒ Yes ⇒ OK button ⇒ Back button (go back one level) ⇒ Diskette symbol to save “empty” message.
- Check dispensing in each program and adjust settings as necessary.
- Perform a function check.
- Back up data via service mode.
Service mode ⇒ Main menu ⇒ Backup ⇒ Backup data for EW ⇒ Overwrite existing data? ⇒ Diskette symbol to overwrite EW data.
- Switch off the machine to exit service mode.

Finish up:

- Check all water connections for leaks.
- Re-install the chassis lid, the lid on the water side, the liquid dispensing cover and the connection box cover.
- Perform an electrical safety check.
- Reconnect the machine to the power supply.

am