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

M.-Nr. 09149681


1 von 26

| | | | | | |
|-----------|---------------------------------------------------------------------------------|-----------|----------------------------------------------------------------|-----------|-------------------------------------------------------------------------------|
| de | Flüssigdosierung, Pumpe DOS-P2, Gehäuse DOS-G2, Ständer DOS-S | en | Liquid dispensing, pump DOS-P2, housing DOS-G2, rack DOS-S | el | Δοσομέτρηση υγρού απορρυπαντικού, αντλία DOS-P2, περίβλημα DOS-G2, βάση DOS-S |
| es | Dosificación de detergente líquido, bomba DOS-P2, carcasa DOS-G2, soporte DOS-S | fr | Dosage liquide, pompe DOS-P2, boîtier DOS-G2, montant DOS-S | it | Dosaggio liquido, pompa DOS-P2, involucro DOS-G2, supporto DOS-S |
| nl | Vloeistofdoserend, pomp DOS-P2, behuizing DOS-G2, standaard DOS-S | pt | Dosagem de líquidos, bomba DOS-P2, caixa DOS-G2, suporte DOS-S | | |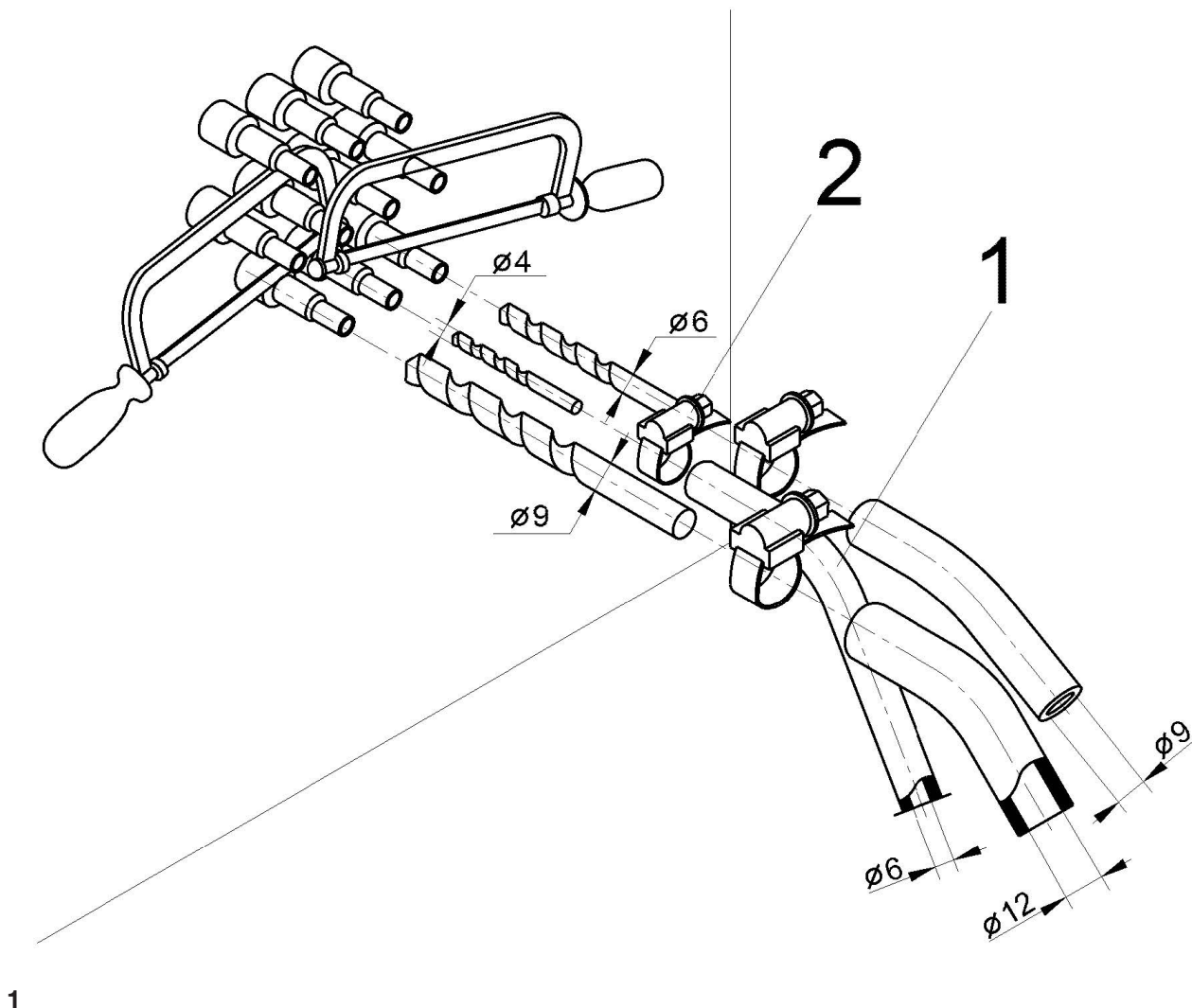


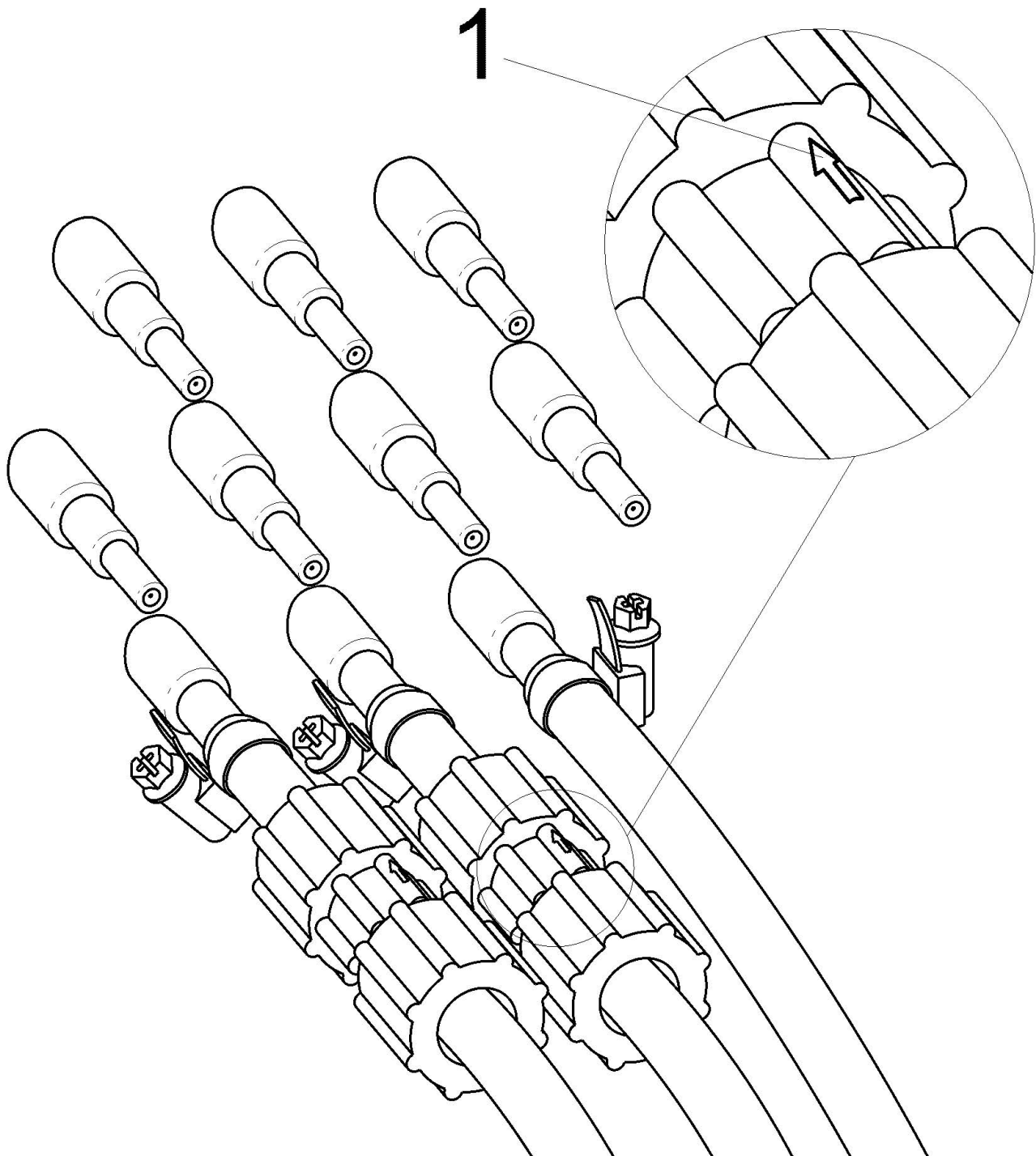
2

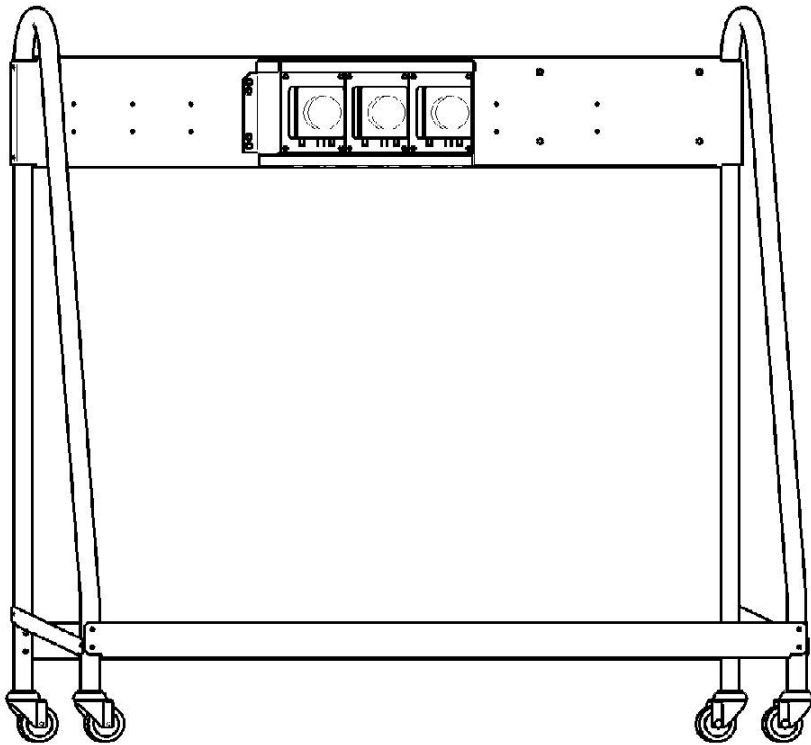3

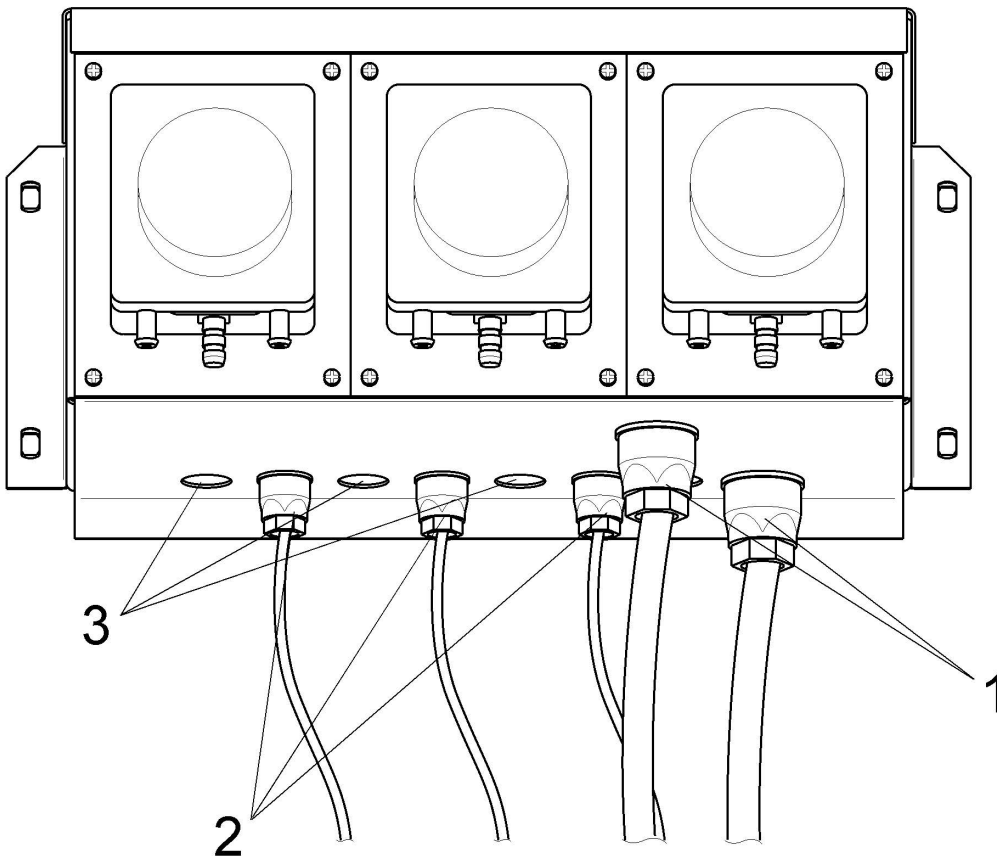4

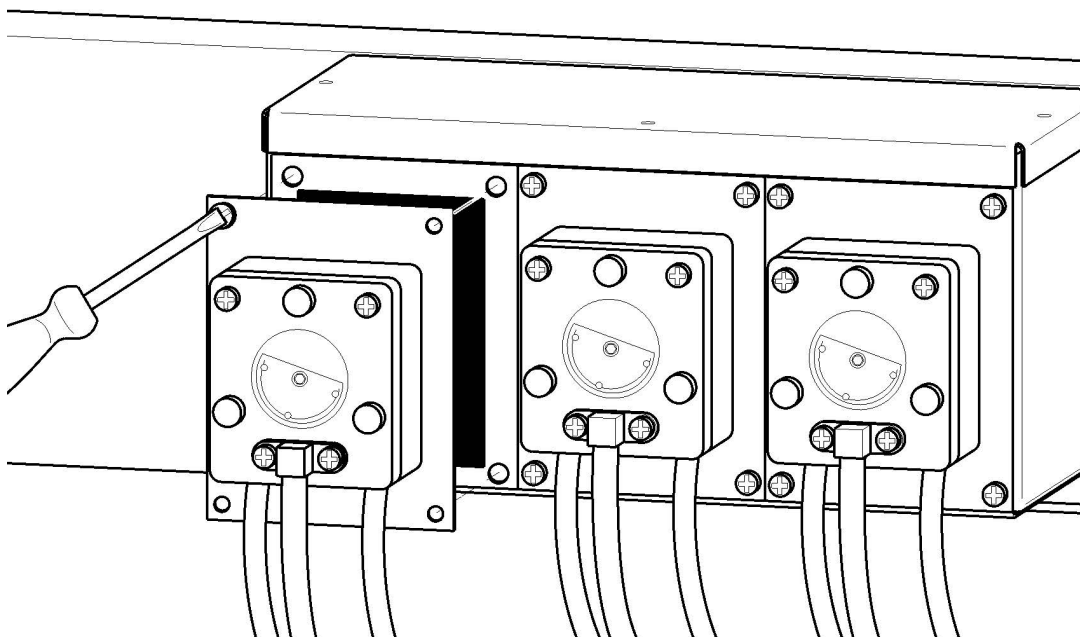5

de

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de

en

Liquid dispensing, pump DOS-P2, housing DOS-G2, rack DOS-S

Reason: these instructions outline the process of installing the dispenser pump DOS-P2, the dispenser pump housing DOS-G2 and the dispenser pump rack DOS-S.

Parts required

| No. | Mat. no. | Designation |
|-----|----------|-----------------------------------------------------------------------|
| | 05573160 | DOS-P2 peristaltic pump |
| | 07569230 | DOS-GP3 Dispenser unit |
| | 04814240 | DOS-S dispensing pump rack for max. 3 pumps DOS-R Non-return valve |

The DOS-P2 dispensing pump fitting kit contains one dispensing pump including accessories for connection to a washing machine:

- Dispensing pump
- Suction lance
- Hoses, clips
- Connection parts for connection to the machine
- Fitting instructions - Pump hose
- Fitting instructions - Liquid dispensing

The DOS-R non-return valve installation kit is required with the use of acidic agents, bleaching agents or similar chemicals, in order to prevent backflow of chemicals (e.g. when containers are changed) into the suction pipework. Non-return valves must not be used with alkaline agents or with conditioning agents or liquid starch.

Check with the detergent supplier.

Contents:

- Non-return valve
- Fitting instructions - Non-return valve

The DOS-G2 dispensing pump housing fitting kit is required for fitting DOS-P2 dispenser pumps. The dispensing pump housing is suitable for a maximum of 3 dispenser pumps.

Contents:

- Housing with lid
- Connection cables
- Wiring material, fastenings
- Wiring diagram - Liquid dispensing, WS 5101 - WS 5323
- Wiring diagram - Liquid dispensing, PW 6101 - PW 6323
- Wiring diagram - Liquid dispensing, PW 6107 - PW 6207
- Wiring diagram - Liquid dispensing, PW 811 - PW 818

The DOS-S dispensing pump stand fitting kit is suitable for the optional securing of the DOS-G2 dispensing pump housing for a maximum of 3 dispenser pumps.

Contents:

- Fastenings for the DOS-G2 dispensing pump housing and the spray proofing pump.

The DOS-GP3 dispenser unit (mat. no.: 07569230) contains the following components:

Included parts

| No. | Mat. no. | Designation |
|-----|----------|--------------------------------------|
| 1 | 07569230 | DOS-G2 pump housing for max. 3 pumps |
| 3 | 05573680 | Dispensing pump |

Parts required for **USA** version:

Parts required

| No. | Mat. no. | Designation |
|-----|----------|------------------------------|
| | 05759720 | DOS-P2 peristaltic pump, USA |
| | 07569230 | DOS-GP3 Dispenser unit |

Parts required for **China** version:

Parts required

| No. | Mat. no. | Designation |
|-----|----------|-----------------------------|
| | 10567550 | DOS-P2 peristaltic pump, CN |
| | 10567650 | DOS-GP3 dispenser unit, CN |

Parts required for **Russia** version:

Parts required

| No. | Mat. no. | Designation |
|-----|----------|-----------------------------|
| | 10864740 | DOS-P2 peristaltic pump, RU |
| | 10864750 | DOS-GP3 dispenser unit, RU |

Depending on the application, one of the following installation kits is also required:

Additional parts required

| No. | Mat. no. | Designation |
|-----|----------|-------------------------------------------------------------------------------------------------------------------------------------|
| | 09079440 | MF01 10-20 multifunction module PW 6xx7 (installation kit – MF01 multifunction module for dispensing pumps 1–6) |
| | 09079450 | MF02 10-20 multifunction module PW 6xx7 (installation kit – MF02 multifunction module for dispensing pumps 7–12) |
| | 09079460 | MF03 10-20 multifunction module PW 6xx7 (installation kit – MF03 multifunction module for dispensing pumps 7–11) |
| | 09079470 | MF05 10-20 multifunction module PW 6xx7 (installation kit – MF05 multifunction module for dispensing pumps 1–3 and water recycling) |

The “Miele Service Documentation (MSD)” service application is required when carrying out maintenance work on the appliance

The safety instructions and warnings in the Miele Service Documentation (MSD) for the corresponding model must be observed.

Work on the appliance must be carried out in accordance with the work instructions.

Installing module for liquid dispensing

- ✚ Remove the machine lid; see the model-specific TSD.
- ✚ For each dispensing pump saw off a connection stub on the mixer box to suit the internal diameter of the hose. When doing so start from underneath. See Fig. 1.
- ✚ Use a 4 mm drill bit to drill a hole in the cut-off stub. See Fig. 1
- ✚ Push hose (1) onto the connection stub, and secure it with a hose clip (2). See Fig. 1.
- ✚ If using a non-return valve cut the hose before the connection stub.
- ✚ Connect both ends of the hose to the non-return valve and tighten the nuts. Observe the direction of flow in the non-return valve. See Fig. 2.
- ✚ Install dispensing pump housing DOS-G2 in its intended location (e.g. the DOS-S dispensing pump rack). See Fig. 3.
- ✚ Fit counter nuts to the cable feed-throughs (1) and (2) in the base of the DOS-G2 dispensing pump housing. See Fig. 4.
- ✚ Close superfluous openings (3) with stoppers. See Fig. 4.
- ✚ Connect the connection cable to the dispensing pump and suction lance according to the wiring diagram.
- ✚ Fit the dispenser pumps DOS-P2 in the dispenser pump housing DOS-G2 and screw them in position. See Fig. 5.
- ✚ Connect the hoses to the dispensing pumps with hose clips.
- ✚ If using less than 3 DOS-P2 dispensing pumps, the vacant pump slot should be closed with the cover provided.
- ✚ Fit the casing lid with 3 self-tapping screws onto the dispensing pump housing DOS-G2.

NOTE

Electromagnetic radiation

Damage to electrical appliances

- ✚ The connection cables for the dispensing pumps and suction lances must be laid separately in order to avoid interference in the control circuit (electromagnetic compatibility). For safety reasons the connection cables must not exceed 4.5 m in length.

- ✚ Pass the connection cables through the cable throughfeeds on the connection strip to the built-in multifunction module and connect them.
- ✚ Refit the machine lid.
- ✚ Connect the machine to the power supply and switch it on.
- ✚ Reconnect the machine to the gas supply if there is one and it has previously been disconnected.
- ✚ Carry out appropriate electrical safety and functional tests.
- ✚ Carry out a functional check only with the dispensing pump system primed and filled.

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