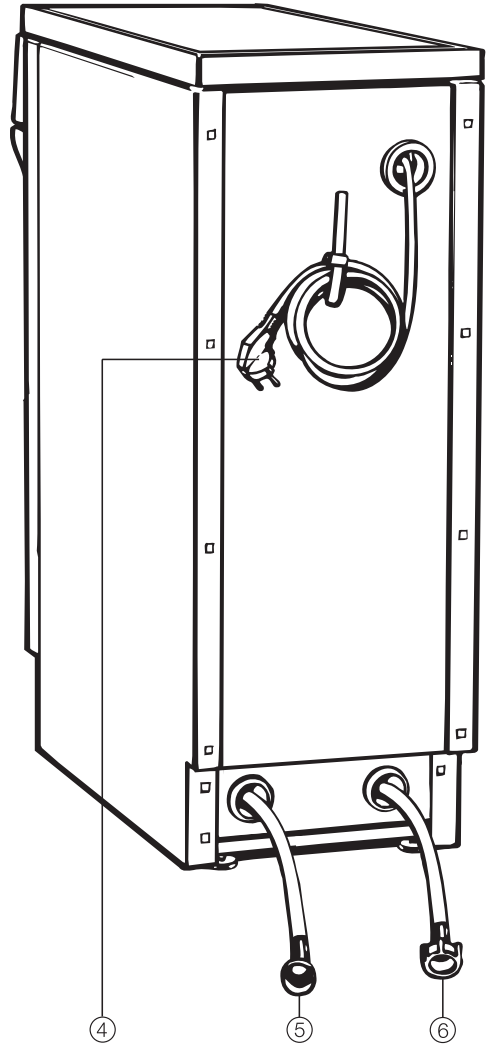
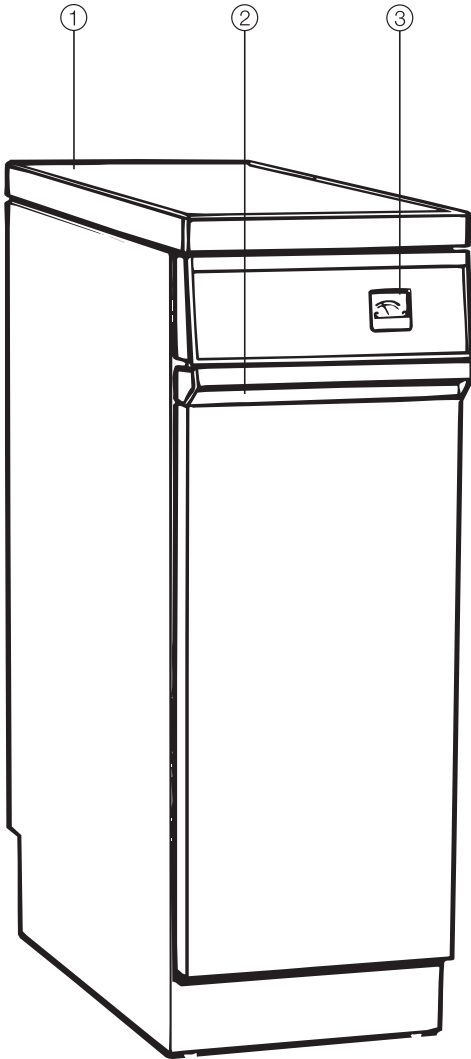
- de** Gebrauchsanweisung Aqua-Purificator
G 7895 und Leitwertmessmodul C
- en** Operating instructions Aqua Purificator
G 7895 and Conductivity Module C
- fr** Mode d'emploi Déminéralisateur G 7895
et module de mesure de conductance C
- it** Istruzioni d'uso Aqua-Purificator G 7895
e Modulo misurazione conducibilità
dell'acqua C
- nl** Gebruiksaanwijzing Aqua-Purificator
G 7895 en module C voor meting
geleidbaarheid (met controlecertificaat)

Inhalt

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
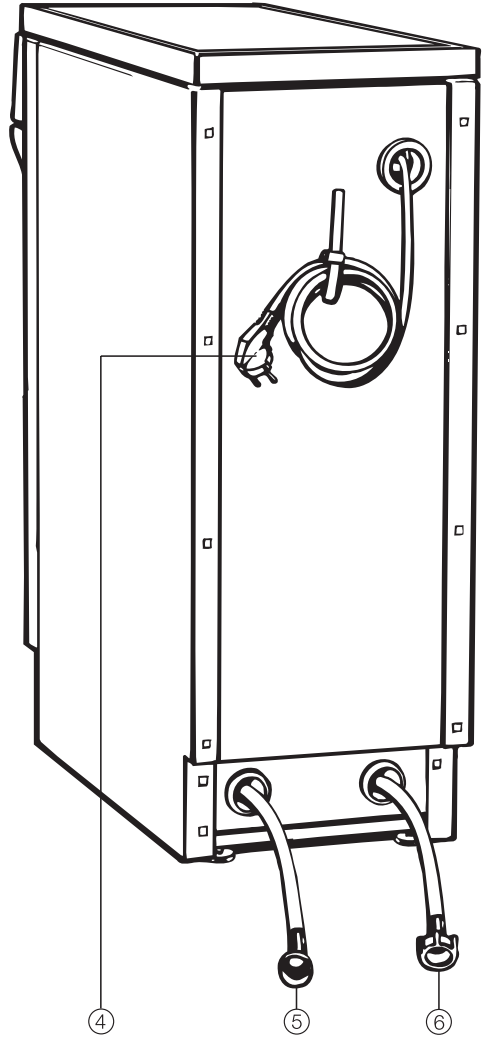
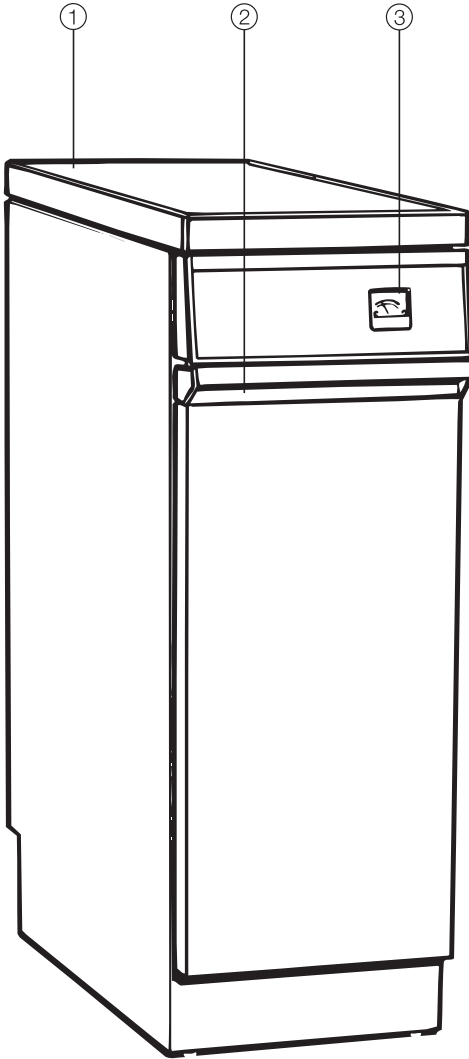
en - Guide to the appliance

- ① Lid
(remove if building under)
- ② Handle for opening drop-down door
- ③ Conductivity display

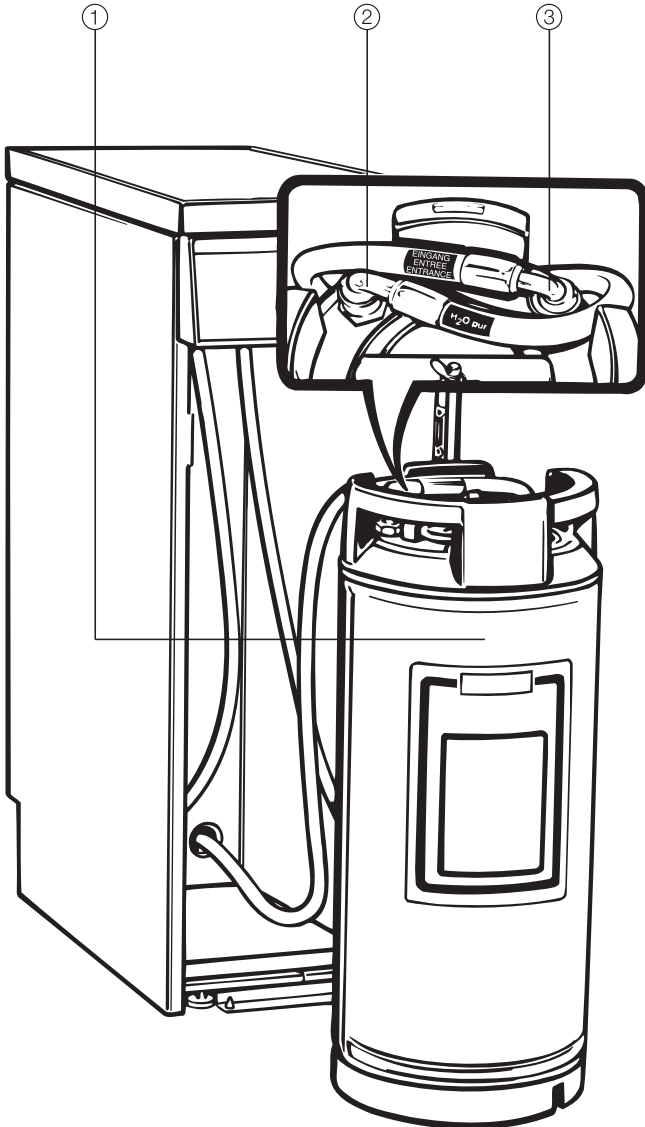- ④ Mains connection cable
- ⑤ Inlet hose (blue)
- ⑥ "H₂O pur" hose for AD water intake
(green)

- ① Water softener canister
- ② Outlet connection
- ③ Inlet connection

en - Warning and Safety instructions

The Aqua Purificator G 7895 and Conductivity module C comply with all relevant safety requirements. Incorrect use can, however, lead to personal injury and damage to property.

To avoid the risk of accidents and damage to these appliances please read these instructions carefully before starting to use them. They contain important notes on the installation, safety, operation and care of the appliances.

Keep these instructions in a safe place for reference, and pass them on to any future user.

Correct application

▶ The Aqua Purificator G 7895 and Conductivity module C are designed for specialised applications only, and must only be used as described in these instructions. Alterations or conversions to the appliances, or using them for purposes other than those for which they are designed are not permitted and could be dangerous.

The manufacturer cannot be held liable for damage caused by improper or incorrect use.

▶ Personnel operating the appliances should be trained regularly. Untrained personnel should not be allowed access to the appliances or their controls.

▶ These appliances are not intended for use by persons (including children) with reduced physical, sensory or mental capabilities, or lack of experience or knowledge, unless they have been shown how to use them correctly by a person responsible for their safety.

Technical safety

▶ The Aqua Purificator G 7895 and Conductivity module C must be checked for any visible damage before they are installed. Do not install and use a damaged appliance. A damaged appliance is dangerous.

▶ The G 7895 should be commissioned and then maintained only by a Miele authorised and trained service technician or other suitably qualified and competent approved person.

▶ The electrical safety of this appliance can only be guaranteed if connected to a correctly installed earthing system on site. It is most important that this basic safety requirement is present and regularly tested, and where there is any doubt the on-site wiring system should be inspected by a qualified electrician. The manufacturer cannot be held liable for damage or injury caused by missing or defective earthing, such as electric shock.

en - Warning and Safety instructions

▶ A damaged appliance is dangerous. Switch off at the mains immediately and call the Miele Service Department or an authorised Miele Service Dealer.

▶ To avoid the risk of damage to the appliance, make sure that the rating on the data plate corresponds to the voltage and frequency of the on-site electrical supply, before connecting it to the mains supply. Consult a qualified electrician if in any doubt.

▶ Installation, maintenance and repairs should only be carried out by a suitably qualified and competent person in accordance with local and national safety regulations. Repairs and other work by unqualified persons could be dangerous and the manufacturer will not be held liable.

Ensure current is not supplied to the appliance until after installation, maintenance or repair work has been carried out.

▶ If the supply cord is damaged, it must be replaced by a Miele approved service technician to avoid a hazard.

▶ Never open the appliance casing. Tampering with electrical connections or components and mechanical parts is highly dangerous to the user and can cause operational faults.

▶ Faulty components must only be replaced by genuine Miele original parts. The manufacturer can only guarantee the safety of the appliance when Miele replacement parts are used.

▶ Do not connect the appliance to the mains electricity supply by a multi-socket unit or an extension lead. These do not guarantee the required safety of the appliance (e.g. danger of overheating).

▶ A damaged appliance is dangerous. Switch off at the mains immediately and call the Miele Service Department or an authorised Miele Service Dealer.

▶ During installation, maintenance and repair work, the appliance must be disconnected from the mains electricity supply. It is only completely isolated from the electricity supply when:

- the mains fuse is disconnected,
- the screw-out fuse is removed (in countries where this is applicable),
- it is switched off at the isolator.
- it is switched off at the wall socket and the plug removed.
- it is switched off at the fused spur connection unit.

Ensure that current is not supplied to the appliance until after maintenance or repair work has been carried out.

▶ In order to fulfill local and national water regulations, a suitable non-return valve must be installed between the mains water connection and the water softener canister.

The non-return valve must be installed at least 200 mm above the highest water level.

en - Warning and Safety instructions

Correct use

▶ Please follow the advice on installation in these instructions and the separate Installation Instructions.

▶ Water in the appliance is not suitable for drinking.

▶ Only tested and approved single use resins may be used for resin exchange.

▶ Never clean the appliance or near vicinity with a water hose or a high pressure hose.

▶ To avoid any corrosion damage, ensure that solutions or steam containing hydrochloric acid do not come into

contact with the steel outer casing of the appliance.

▶ While the appliance is under guarantee, repairs should only be undertaken by a service technician authorised by the manufacturer. Otherwise the guarantee will be invalidated.

▶ This appliance must only be used in conjunction with machines specified by Miele as being suitable for such usage.

Installation

▶ Do not install the Aqua Puricator in an area where there is any danger of explosion or of freezing conditions.

▶ This equipment may only be used in mobile installations such as ships if a risk assessment of the installation has been carried out by a suitably qualified engineer.

Safety with children

▶ To avoid the risk of injury, keep children away from this equipment at all times. To avoid the risk of injury, do not allow them to play with it or use the controls. They will not understand the potential dangers posed by it. There is a danger that they could injure themselves or get into the appliance and lock themselves in.

▶ Older children may only use this equipment if its operation has been clearly explained to them and they are able to use it safely, recognising the dangers of misuse. They must be aware of the potential dangers caused by incorrect operation.

Disposing of old appliances

► For environmental and safety reasons ensure the appliance is completely drained of any residual water and resin. Observe safety regulations whilst doing so. Make the door lock inoperative, so that children cannot accidentally shut themselves in. Then make appropriate arrangements for the safe disposal of the appliance.

► Single use resin and resin that cannot be reactivated can be disposed of with the normal waste disposal.

The manufacturer cannot be held liable for damage caused by non-compliance with these Warning and Safety instructions.

en - Caring for the environment

Disposal of the packing material

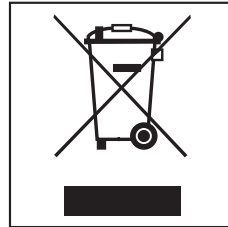
The packaging is designed to protect the appliance during transportation. It has been selected from materials that are environmentally friendly for disposal and that can be recycled.

Recycling saves on the use of raw materials and reduces the amount of waste going to landfill.

Ensure that any plastic wrappings, bags etc. are kept out of the reach of babies and young children. Danger of suffocation. Depending on country your dealer may also be able take packaging back from you.

Disposal of your old appliance

Electrical and electronic appliances often contain materials which, if handled or disposed of incorrectly could be potentially hazardous to human health and to the environment. They are, however, essential for the correct functioning of your appliance. Please do not therefore dispose of your old appliance with your general waste.

Please dispose of it at your local community waste collection/recycling centre, or contact your dealer for advice.

Please store it out of reach of children whilst awaiting disposal.

en - Description of the appliance

The **Aqua Purificator G 7895** can be fitted with one of two different water softening canisters:

– Water softener canister E 310

E 310 canisters are filled with reusable resins, which can be regenerated once depleted. The whole canister must be sent to a regeneration station for reprocessing.

For the address, contact your Miele dealer.

– Water softener canister E 318

E 318 canisters are supplied empty, and have to be filled on site with 20 l single-use resins (see "Filling an E 318 canister with single-use resin").

Once depleted, these resins have to be replaced by new single-use resins.

Single use resin and resin that cannot be reactivated can be disposed of with the normal waste disposal.

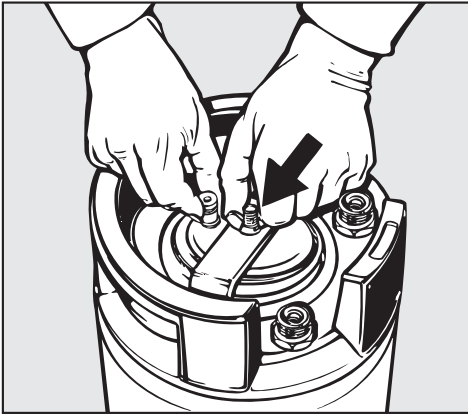
To easily fill and re-fill an E 318 canister, we recommend using the E 316 Miele Resin Exchange Set consisting of:

- A 30 l plastic barrel with lid and funnel.
- 2 x E 315 bags of single-use resin (each 10 l), with a filter bag for the spent resin.


Additional 20 l bags of single-use resin can also be ordered as a set under code E 315.

en - Fill the E 318 canister with E 315 single-use resin.

Open the lid

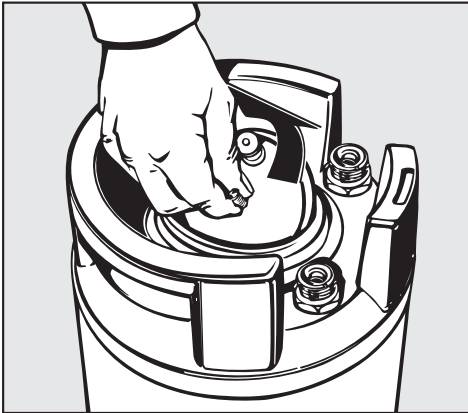
- Undo the wing nuts, holding the lid to prevent it from falling into the canister.
- Remove the clamp.

- Lift the lid.

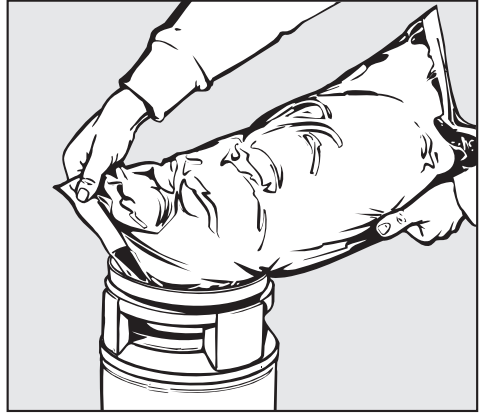
Fill with resin

- Place the funnel in the canister.
- Cut a corner from the top of each bag of resin, and empty them into the canister.
- Wipe up any resin from around the opening.
- Replace the lid, holding the vent firmly and secure with the clamp and wing nut.

en - Connecting the E 310 or E 318 canister to the unit.

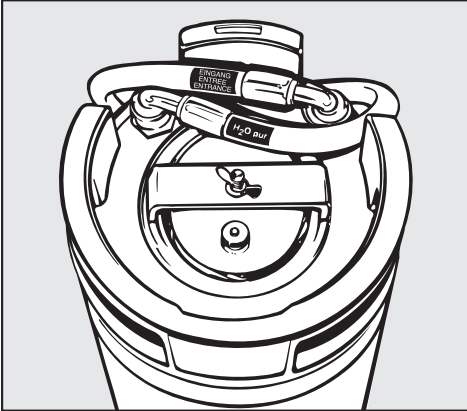- Open the front of the Aqua Purificator.
- Unscrew the safety caps from the connection points on the canister.

- Attach the blue inlet hose to the "inlet", and the green H₂O pur hose to the "outlet".
- Arrange the hoses as shown in the diagram.

en - Connecting to the Aqua Purificator

- Connect the green hose marked "H₂O pur" to the AD water solenoid valve or to the AD inlet connection hose on the machine.
- Connect the blue "Inlet" hose to the on-site cold water stopcock.

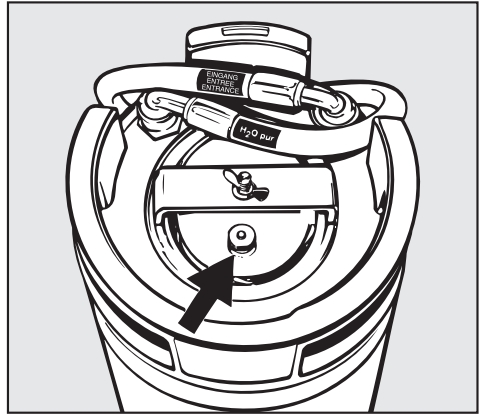
See enclosed installation instructions for further details.

Important

The water pressure (flow rate) must be between 250 and 1000 kPa (2.5 -1 10 bar).
The water temperature must not exceed 20 °C.

- The plug must be correctly inserted in the socket and switched on for the conductivity display to work.

Vent the canister

- Open the vent (see arrow) by turning it anti-clockwise.
- Open the cold water stopcock slowly until water starts to come out of the vent.
- Then close the vent.

Install the canister

- Fit the connected canister in position at **the back** of the Aqua Purificator. The area at **the front** can be used to store a spare canister.
- The Aqua Purificator is ready for use once the shut-off valve is open.

The conductivity display shows the conductivity of the fully demineralised water, which is measured in $\mu\text{S}/\text{cm}$ (microsiemens per cm). This enables the softened water to be quality controlled.

If the minimum water quality required is not being achieved, the canister will need to be reactivated.

- **Canister E 310:** Canisters filled with reusable resins must be sent to a regeneration station for reactivation.
- **Canister E 318:** Replace the resin. Instructions on how to do this are supplied with the E 316 Resin Exchange Set.

Removing the E 310 or E 318 canister

- Turn off the cold water stopcock.
- Open the stopcock for manual AD water (if present), or
- Open the vent on the canister to release the pressure.
Mop up any spilt water with a cloth.
- On Miele washer-disinfectors equipped with **programme card controls** insert programme card "Abpumpen/ETT" (Drain/ETT) to pos. 0 to release the build-up of pressure, and then switch the machine on for a few seconds.
- Disconnect the hoses.
- Up-end the canister and empty out any residual water.

Sending off the E 310 canister (with reusable resin)

- Screw the plastic safety caps to the "inlet" and "outlet" connection points on the canister before sending it away for regeneration.

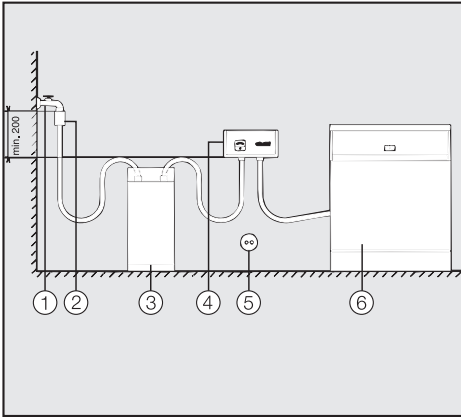
To avoid the risk of frost damage

Fill the canister with approx. 500 g cooking salt before sending it off. Secure the lid with wire.

en - Conductivity module

The conductivity module fulfills the same role as the integrated conductivity display on the Aqua Purificator G 7895.

Measuring the conductivity level of the water enables the fully demineralised water to be quality controlled.

- The conductivity module is always connected behind the water softener canister in the following order:
Stopcock ① - Non-return valve with air gap ② - Water softener canister ③ - Conductivity module ④ - Machine ⑥.
- It can be installed on a wall or in a base unit (it must **not** be installed on the floor or under taps).
- Electrical connection ⑤ via plug and fused socket.

See enclosed installation instructions for further details.

 All electrical work must be carried out by a suitably qualified and competent person in accordance with current local and national safety regulations (BS 7671 in the UK).

- Ensure power is not supplied to this appliance until after it has been installed.
- Connection should be made via a suitable switched socket which is easily accessible for servicing and maintenance work after the appliance has been installed. An electrical safety test must be carried out after installation and after any servicing work. When switched off there must be an all-pole contact gap of at least 3 mm in the mains connection switch.
- The mains connection cable may only be replaced by an original Miele spare part or an appropriate cable with **wire end ferrules**.
- For technical data, see data plate.

The machine must only be operated with the voltage, frequency and fusing shown on the **data plate**.

Non-rewireable plugs BS 1363

The fuse cover must be refitted when changing the fuse, and if the fuse cover is lost, the plug must not be used until a suitable replacement is obtained. The colour of the correct replacement cover is that of the coloured insert in the base of the plug, or the colour that is embossed in words on the base of the plug (as applicable to the design of plug fitted).

Replacement fuses should be ASTA approved to BS 1362 and have the correct rating. Replacement fuses and fuse covers may be purchased from your local electrical supplier.

WARNING: THIS APPLIANCE MUST BE EARTHED

en - Certification

The Miele Aqua Purificator G 7895 with water softener canister E 310 or E 318 must be safety tested after commissioning and after any maintenance work in accordance with local and national safety regulations on the use of pressurised containers.

– A - Check that the following are all present

At the installation site, make sure the following are all present (see sticker on canister):

1. Certification of the pressure test by the manufacturer of the pressurised canister.
2. Instructions on how to use the pressurised canister.
3. Technical data for the pressurised canister.

– B - Equipment check

1. Make sure all components are present:
 - Pressure hose from supply to canister
 - Pressure hose from canister to conductivity sensor
 - Pressure hose from conductivity sensor to machine **or**
 - Pressure hose from conductivity sensor to machine supplying the water
 - Water supply (on request)
 - Cartridge
2. Check everything works correctly

– C - Test of installation requirements

1. The pressurised canister is housed in the Miele Aqua Purificator G 7895.
2. The Miele Aqua Purificator G 7895 is designed so that it is accessible from the front.

– D - Regular testing according

Points A, B and C must be checked by a Miele Service Technician every 5 years.

Miele Service Department

Date _____

(Signature)

Should you have any difficulties please contact your

- Your Miele dealer, or
- the Miele Service Department
(see back cover for contact details).
- Please quote the model and serial number of your appliance.

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